

**WHITE LABS® ANALYTICAL SERVICES**

# **BIG QC DAY**

## **- TEST KIT -**

**STOP  
BEFORE  
PROCEEDING**

1. Please read all instructions before starting.
2. Register samples following instructions.  
Samples will not be processed if not registered.
3. Return samples by August 23, 2019, to get results completed by September 6, 2019.

### **YOUR KIT INCLUDES THE FOLLOWING ITEMS FOR COLLECTION AND SHIPMENT OF TWO BEER SAMPLES.**

- (2) 250mL sterile sample containers with unique barcodes
- Return shipping label (except international customers)



**PURE YEAST & FERMENTATION**

# 1

## REGISTER SAMPLES ONLINE

- Log in to your White Labs Sample Analysis Customer Portal at <https://whitelabs.qbench.net>.
- If you did not receive an email invitation to create a login or have forgotten your portal password, please email [analyticalab@whitelabs.com](mailto:analyticalab@whitelabs.com) or call (858) 693-3441 ext. 7387.
- Once you are signed in, click “Register Big QC Day Samples Here” at the top of the page.
- Select the Order ID.

**Order Requests**  
White Labs Tasting Room / Order Requests / Filtered

Show 10 entries

ID	Request Status	Date Received
26405	PENDING CHANGES	07/17/2019 10:30 AM

Showing 1 to 1 of 1 entries

- Insert sample name in the description field for each Sample ID.

Samples

Sample ID	Sample Type	Description	Comments
39141	Beer		+ Add
39142	Beer		+ Add

[Save Samples](#)

- Click “Save Samples.”
- Click “Submit Sample Information” at the top right.

[Submit Sample Information](#)

## 2

## OBTAIN CLEAN SAMPLES

*All samples must be shipped in the supplied numbered containers.*

### **Obtaining a beer sample from a beer bottle:**

- Have barcoded sample container nearby with lid loosened.
- Use isopropanol spray to clean the outside top of beer bottle and bottle opener.
- Open the beer and carefully pour into the sample container. Fill to the 200mL mark.
- Close lid securely.

### **Obtaining a beer sample from tank:**

- Have a barcoded sample container nearby with lid loosened.
- Clean the sample port with a cotton swab soaked with isopropanol.
- Use isopropanol spray to clean outside of the sample port. Flame faucet with portable flame if possible.
- Let approximately 8oz of beer flow through tap before collecting it into the sample container. Fill to the 200mL mark.
- Close lid securely.
- Repeat cleaning procedure once finished taking the sample.

### 3 SEND SAMPLES TO WHITE LABS

- Put containers in the original shipping envelope and tape. Put the return shipping label on shipping envelope.
- Call FedEx for a pick-up or drop off at your nearest FedEx location.

### 4 GET YOUR RESULTS

There are dates throughout the year that we run this test. Please reference our website for current dates at [whitelabs.com/bigqcday](http://whitelabs.com/bigqcday). Samples must be postmarked by the deadline in order to receive results by the specific date.

Once testing is complete, you will receive an email with the reports. Your results will also be available to view via the White Labs Sample Analysis Customer Portal at [whitelabs.qbench.net](http://whitelabs.qbench.net)



[WHITELABS.QBENCH.NET](http://whitelabs.qbench.net)

[WHITELABS.COM](http://whitelabs.com) | [YEASTMAN.COM](http://yeastman.com)